


### Exploded View for 2090

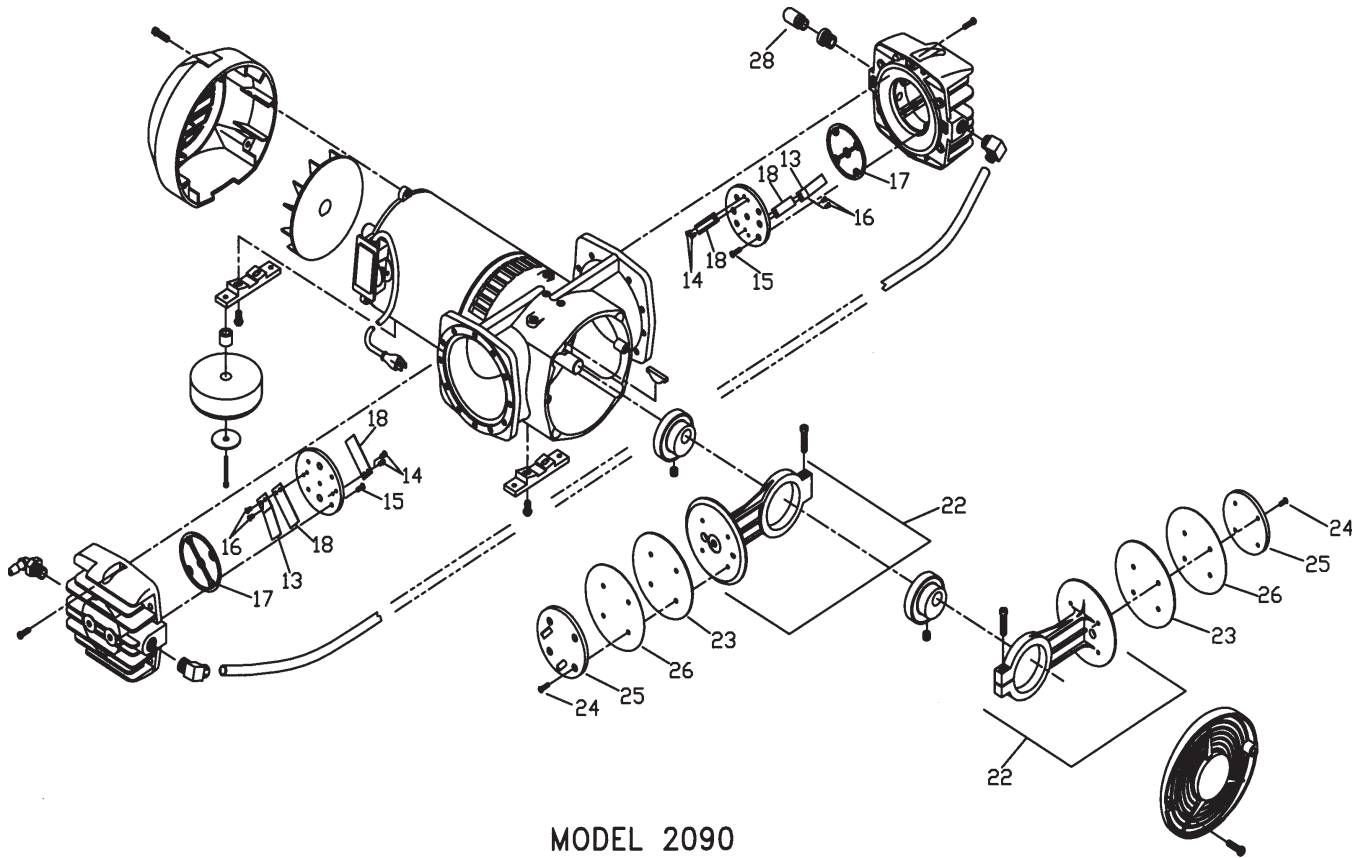

MODEL 2090

67-1341  
SHEET 2 OF 2  
REV. 1 08/01

### Parts List for 2090

Item No.	QTY	Part No.	Description	Diaphragm Service Kit 2090K-01	Valve Plate Service Kit 2090K-01
13	2	See Note 1	Valve Keeper	-	1
14	4	See Note 1	Screw-Valve Flapper-Intake	-	1
15	10	See Note 1	Screw-Valve Plate	-	5
16	4	See Note 1	Screw-Valve Flapper-Exhaust	-	5
17	2	See Note 1	Valve Plate Gasket	-	1
18	4	See Note 1	Valve Flapper-Intake & Exhaust	-	2
22	2	2090K-03	Connecting Rod	-	-
23	2	See Note 1	Diaphragm	1	-
24	8	See Note 1	Screw-Hold Down Plate	4	-
25	2	See Note 1	Hold Down Plate	1	-
26	2	See Note 1	PTFE Liner	1	-
28	1	See Note 1	Silencer	-	-

Note 1: These parts are not available separately, but are supplied within kits.

67-1341  
Sheet 1 of 2  
Rev. 1 08/01