


## Section 10: EXPLODED VIEWS AND PARTS LIST

### 10.10 MODEL 2546

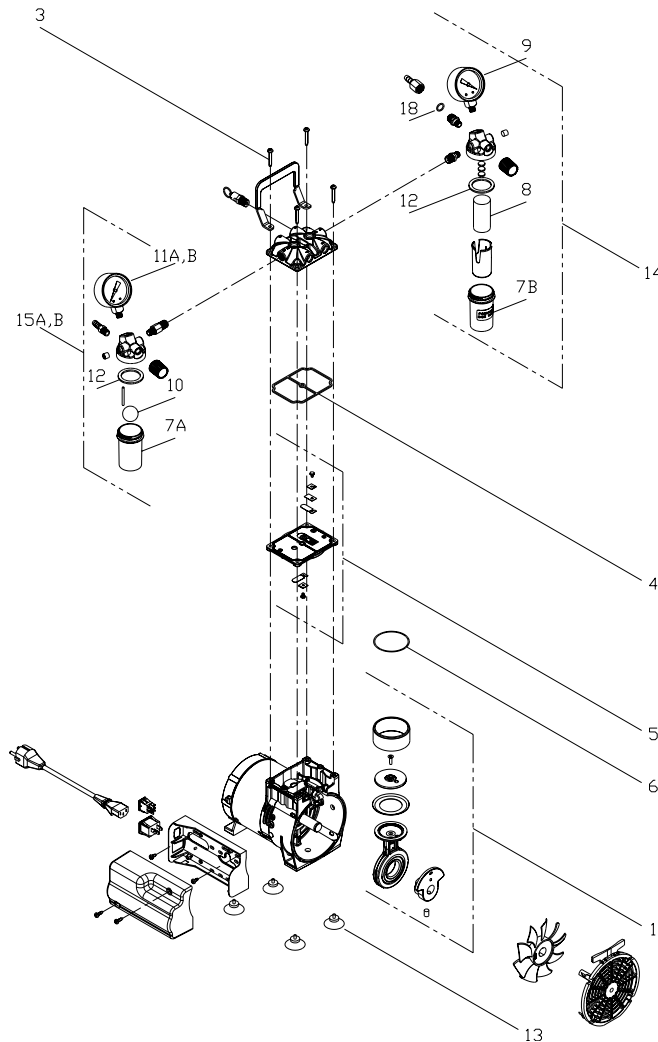

Item Number	QTY	Part Number	Description	Service Kit 2546K-03	Seal Service Kit 2546K-04	Feet Service Kit 2500K-04
1	1	2546K-01	Connecting Rod Assembly	1		
3	4	Note 1	Head Screw	4		
4	1	Note 1	O-Ring Head	1	1	
5	1	2567K-02	Valve Plate Assembly	1		
6	1	Note 1	O-Ring Cylinder	1	1	
7	2	Note 1	Jar Plastic	1	1	
8	1	Note 1	Filter Element	1	1	
9	1	726025	Pressure Gauge			
10	1	Note 1	Ball	1	1	
11A	1	726021	Vacuum Gauge 2546B-01			
11B	1	726024	Vacuum Gauge 2546C-02			
12	2	Note 1	Gasket	1	1	
13	4	Note 1	Rubber Suction Feet			4
14	1	2522K-06	Pressure Regulator Assembly			
15A	1	2522K-05	Vacuum Regulator Assembly 2546B-01			
15B	1	2522K-07	Vacuum Regulator Assembly 2546C-02			
16	1	Note 1	Piston Cup Seal		1	
17	1	Note 1	Retainer Screw		1	
18	1	Note 1	O-Ring	1	1	