


6.14

Parts List and Exploded View for 2561

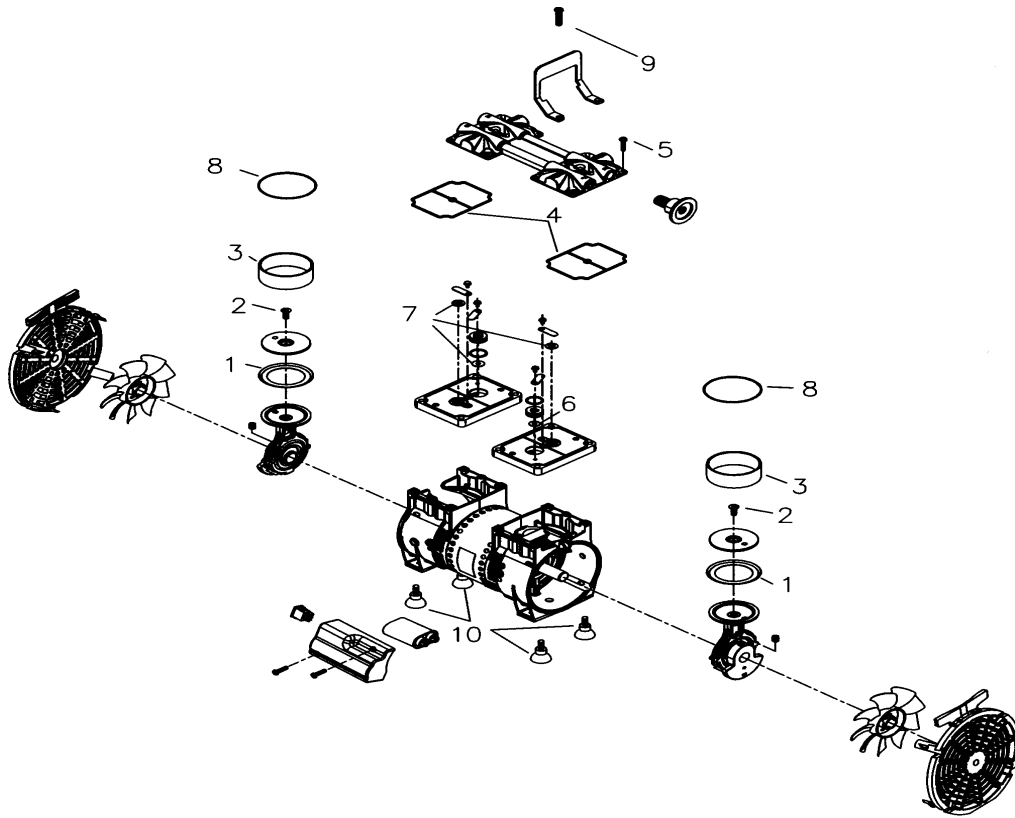

Model 2561

Item No.	Part No.	Description	Qty	Service Kit 2561K-03	Seal Service Kit 2561K-04	Feet Service Kit 2500K-04
1	61-3603 *	Piston Cup	2	2	---	---
2	61-3604 *	Cup Retainer Screw	2	2	---	---
3	61-3441 *	Cylinder Sleeve	2	2	---	---
4	62-1052 *	O-Ring, Gasket	2	2	2	---
5	61-3607 *	Head Screw	6	6	---	---
6	66-0288 *	Intake Valve, Intake Stage	1	1	1	---
7	66-0289 *	Intake/Exhaust Valve, Intake and Exhaust Stage	3	3	3	---
8	62-1050 *	O-Ring, Cylinder Sleeve	2	2	2	---
10	66-0160 *	Feet, Suction Cup	4	---	---	4

* Note: These parts are not available separately, but are supplied within kits shown in last three columns.