


## 6.15 Parts List and Exploded View for Pump Model 2565

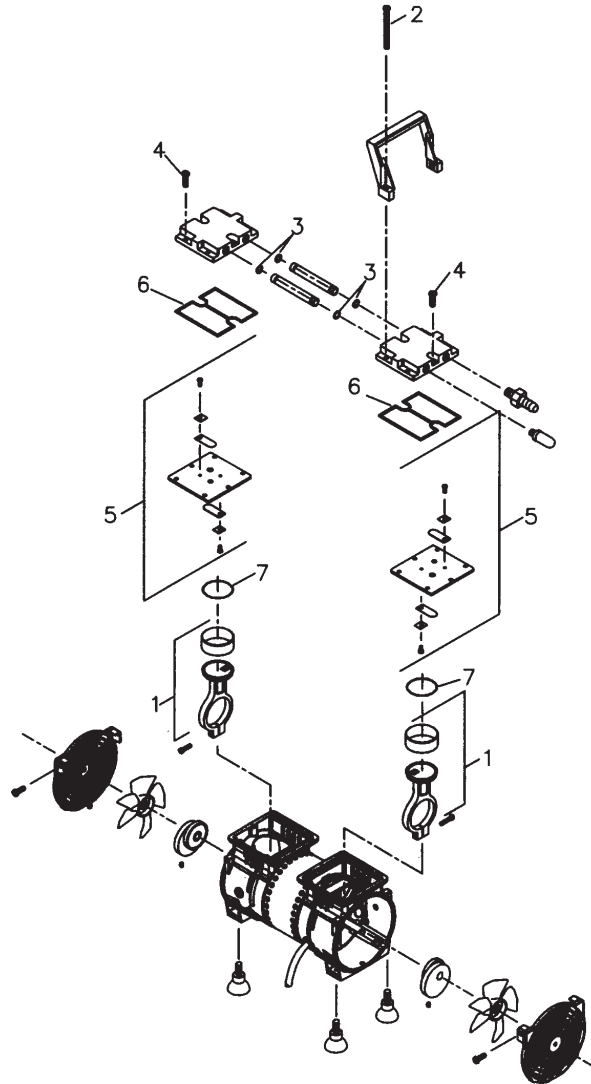

Model 2565

Item Number	Qty.	Part No.	Description	Service Kit 2565K-03	Seal Kit 2565K-04	Feet Kit 2500K-04
1	1	2565K-01	Connecting Rod Ass'y	2	-	-
2	2	Note 1	Handle Screw	2	-	-
3	2	Note 1	O-Ring, Connector Tube	4	4	-
4	10	Note 1	Head Screw	10	-	-
5	2	2565K-02	Valve Plate Ass'y	2	-	-
6	2	Note 1	O-Ring Valve Plate	2	2	-
7	2	Note 1	O-Ring Cylinder	2	2	-
8	1	1412C	Muffler	1	-	-
9	4	Note 1	Rubber Suction Feet	-	-	4

Note 1: These parts are not available separately, but are supplied within the kits shown in last three columns.