


Parts List for Model 8140

Item No.	Qty.	Part Number	Description
1	1	71-1266	Ring for Model 8140
2	1	8120K-01	Vane Kit; 4 Vanes, 1 Gasket
3	1	71-1257	End Plate
4	1	71-1258	Sound Chamber
5	1	71-1259	Gasket
6	1	71-1260	Shroud
7	2	71-1261	Filter Kit; 4 Felts, 1 Ring, 1 Holder
8*	1	71-1268	Exhaust Assembly
9	1	71-1267	Intake Assembly
10	2	71-1275	Filter Jar
11	2	71-1270	Filter Element
12	2	71-1269	Filter Jar Gasket
13	2	62-0337	Brass Nipple
14	1	71-1357	Carbon Wear Plate
15	1	71-1358	Felt Washer

*Model 8140B-15 Does not come with exhaust assembly since it is configured for compressor operation

Exploded View of Model 8140

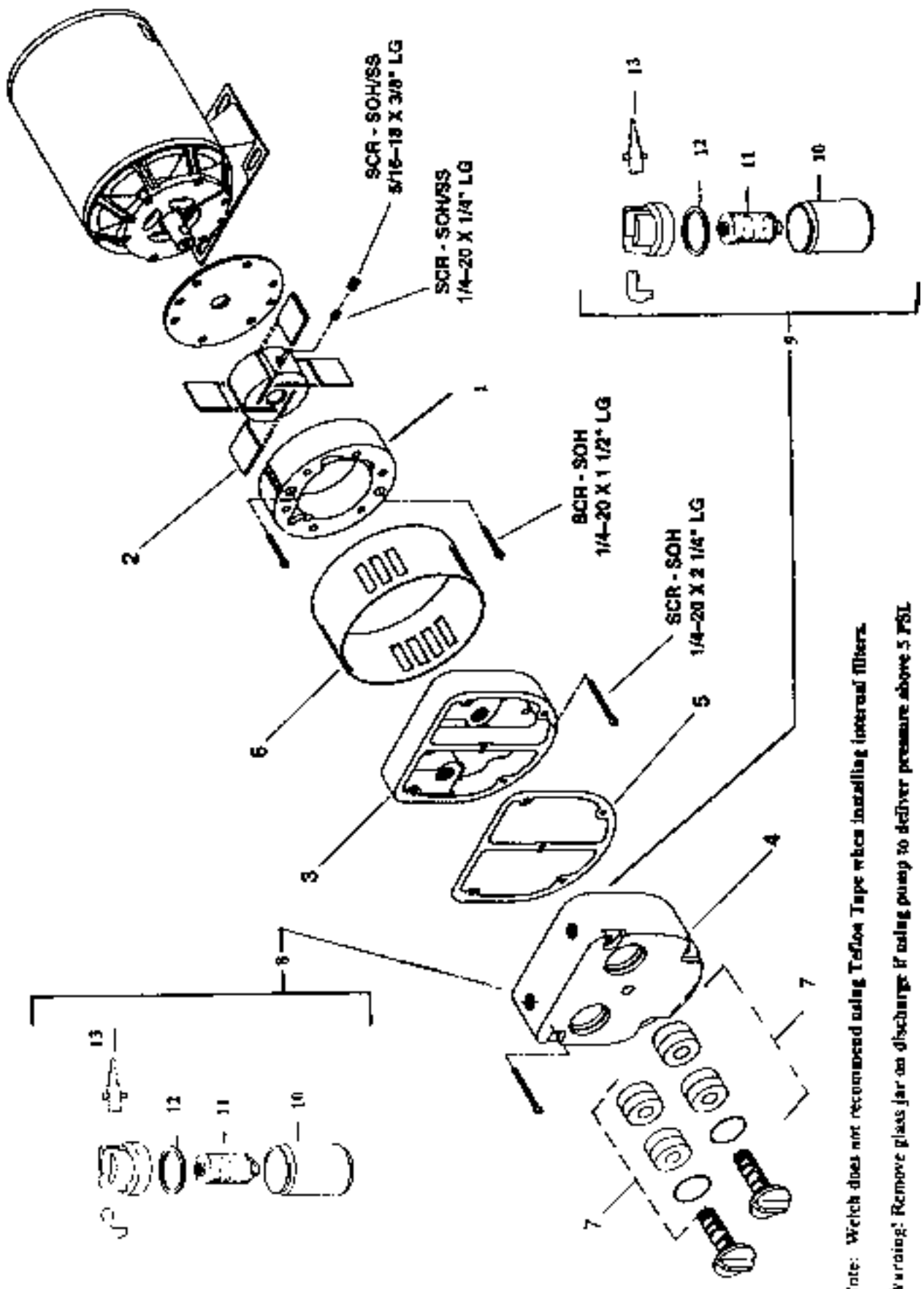

Note: Welch does not recommend using Teflon Tape when installing internal filters.

Warning: Remove glass jar on discharge if using pump to deliver pressure above 5 PSI.